


### ① Quick Coomassie Stain

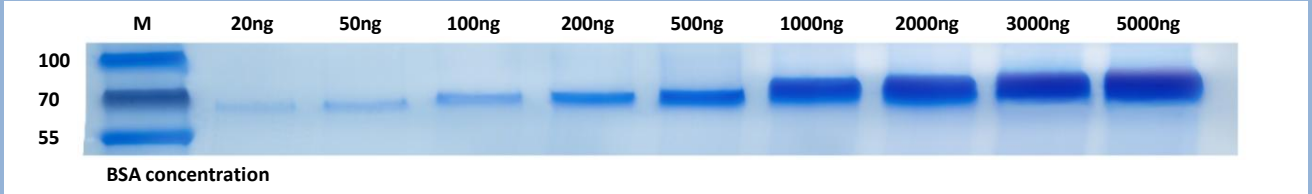


Optimized Post-Stain for complex mixtures, lysates, serum, pure and partially pure proteins

Speed up your SDS PAGE protein staining with:

- FAST • SENSITIVE • SAFE
- 50 times more sensitivity than other stains
- Protein bands down to 5ng
- Rapid 1-step procedure
- Ultra low background

②

**The Only Universal Protein  
Pre-stain on the Market**

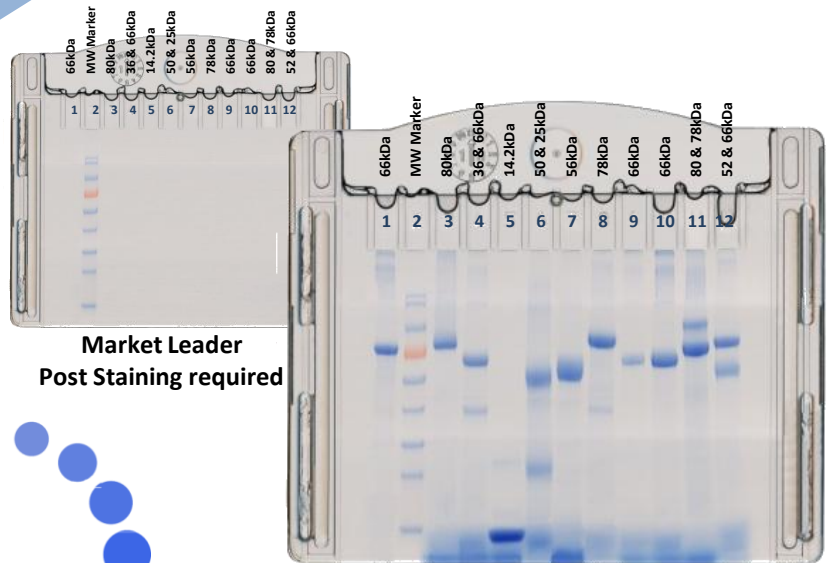

**Realtime**  
stain

Patent Pending

**Optimized for pure and  
partially pure proteins**


Order codes:	Description:	No of gels (10 lanes)
GEN-QC-Stain-1L	Quick Coomassie Stain (1L)	40
GEN-QC-Stain-3L	Quick Coomassie Stain (3L)	120

## HIGH RESOLUTION

Sharp protein bands that you would expect with traditional Coomassie staining

## SENSITIVE

50 x more sensitive than other rapid stains

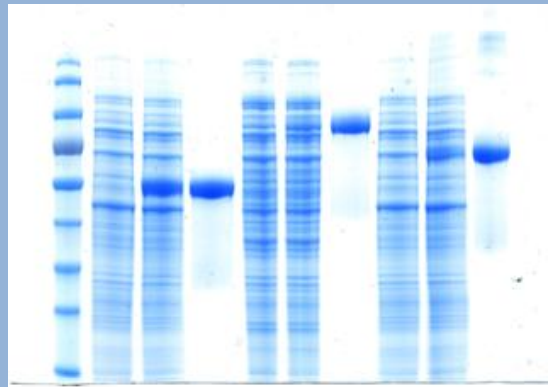

Samples prepared using Tris-Tricine sample loading buffer and run on a 12% NuSep Tris-Glycine SDS PAGE gel. Stained using Quick Coomassie stain.

## RAPID

15 min 1-step stain

## LINEAR RANGE

Very low background to accurately quantify proteins

## DURABLE

Re-usable up to 3 times!  
Also MS compatible

Produced in association with Geron Limited

**Generon**

Dual Trial Pack

## Realtime

View protein band separation in **Realtime** during electrophoresis

Quick Coomassie Stain (30ml)  
Realtime Stain (0.2ml)

## Instant


NO Coomassie post staining required

## Customised

Make in-house protein ladders with your own well-characterised proteins

GEN-QC-RT-Trial

M 6000ng 5000ng 4000ng 2000ng 1000ng 500ng 250ng 100ng 50ng

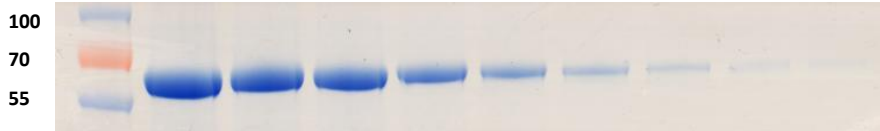

Sensitivity down to 50ng

## UNIQUE

Behaves like Coomassie with NO background

Order codes:	Description:	No. of gels (10 lanes)	No. of lanes
GEN-RT-Stain-2000	Realtime Stain (2ml)	40	400
GEN-RT-Stain-200	Realtime Stain (0.2ml)	4	40